


Unison UK-Phenyl Cadenza CX-C18  
 Unison UK-C8 Cadenza CW-C18  
 Unison UK-C18 Cadenza CD-C18

100 x 4.6 mm


HPLC COLUMN APPLICATION

## Bis [2- [ (oxo) diphenylphosphino ] phenyl ] ether (organic light-emitting diode material)

ビス [ [ 2 - [ (オキソ) ジフェニルホスフィノ ] フェニル ] エーテル  
 (有機エレクトロルミネッセンス材料)

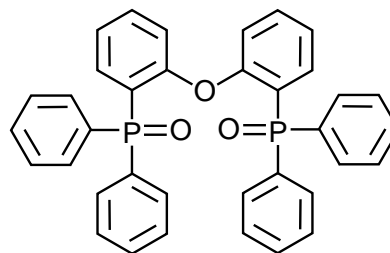

bis [2- [ (oxo) diphenylphosphino ] phenyl ] ether

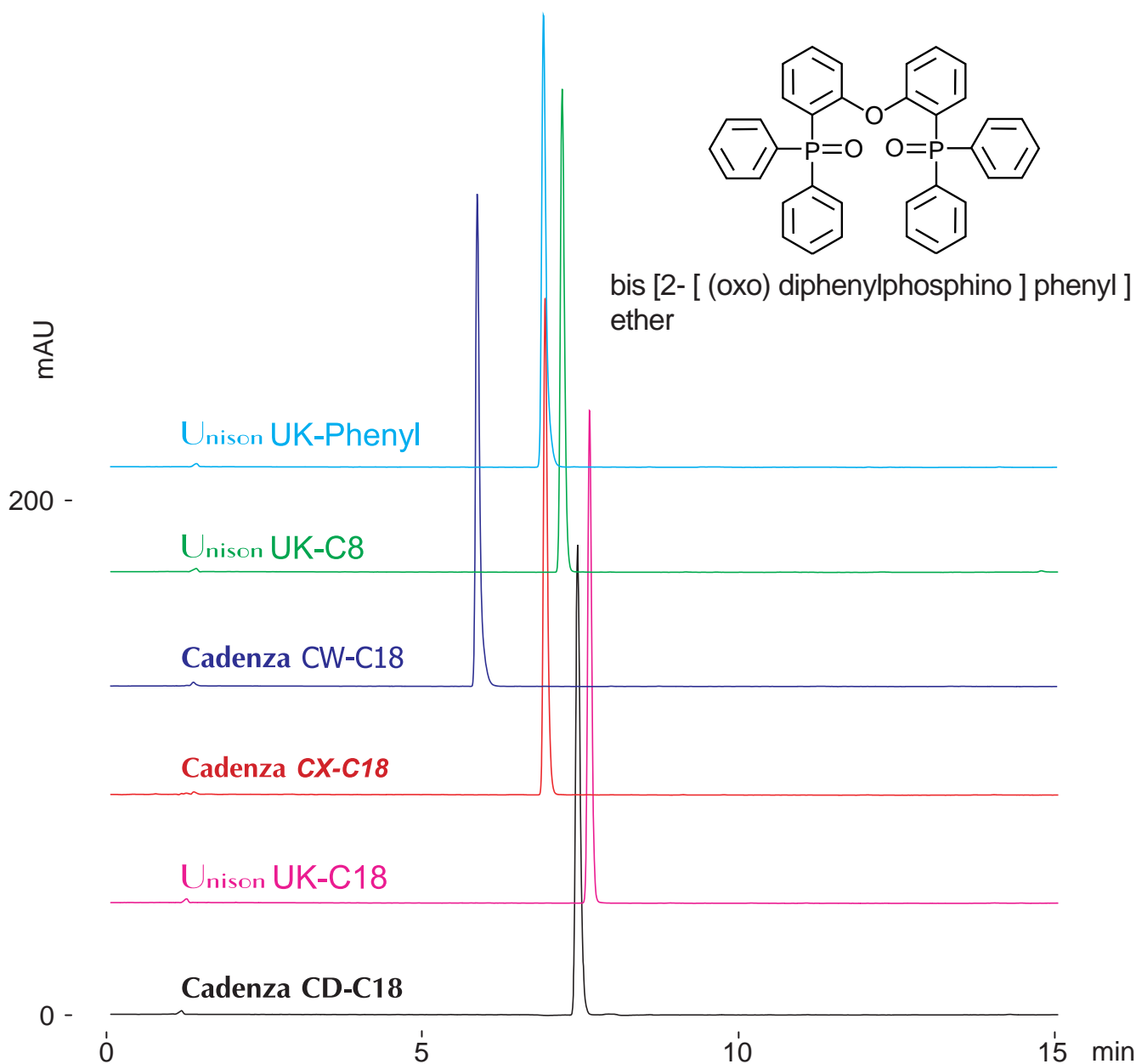

100 x 4.6 mm

A: water / acetonitrile / acetic acid = 90 / 10 / 0.1

B: water / acetonitrile / acetic acid = 10 / 90 / 0.1

25-100%B (0-15min)

1.0mL/min (10MPa), 37°C, 290nm

10µL (0.5mg/mL, water / acetonitrile / acetic acid = 30 / 70 / 1)